METRIA pH tester – pH 600

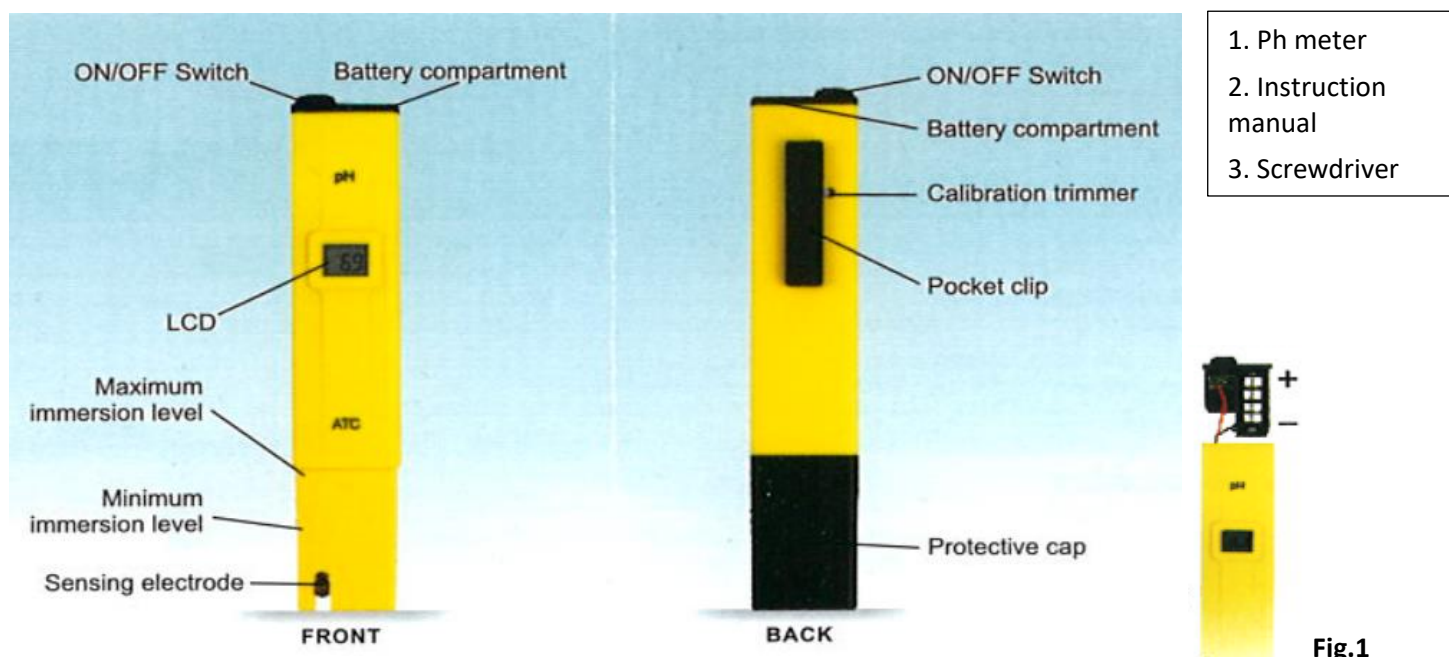
Fig.1

METRIA

pH tester

1. Contains:

pH meter, instruction manual, screwdriver.

2. Operation:

Remove the protective cap

-Do not be alarmed if white crystals appear around the cap. This is normal with pH electrodes and they dissolve when rinsed in water.

-Turn the meter on sliding the switch on the top.

-Stir gently and wait until the display stabilizes. Readings are automatically compensated for temperature variations.

-After use, switch off the pH meter, clean the electrode and put the cap back on.

NOTE: Do not use distilled or deionized water for storage purposes.

Large fluctuations in readings could be due to a lack calibration, electrode dryness or low battery.

3. Calibration:

-Sink the tester up to the maximum immersion level in pH 6.86 buffer solution.

-Allow the readings to stabilize and use the included screwdriver to adjust the calibration nut at the back of the device to read 6.9.

NOTE: Always use fresh buffers for calibration purposes.

4. Battery replacement (Fig.1)

When the meter cannot be switched on or the display fades, open the battery compartment cover and replace all four 1.5V batteries, taking into account the polarity.

Batteries should only be replaced in a safe area using the battery type specified in this instruction manual.

5. Technical data

Resolution	0.1 pH
Accuracy	± 0.1 pH
ATC	0 to 50° C (32 to 122° F)
pH scale:	0.0 to 14.0
Environment	0 to 50° C (32 to 122° F) RH 95% max
Battery type	4 x 1,5 V alkaline batteries
Battery life	Approx. 150 h of use
Dimensions	152 x 30 x 32 mm (5.9x1.2x0.8 ")
Weight	50 g

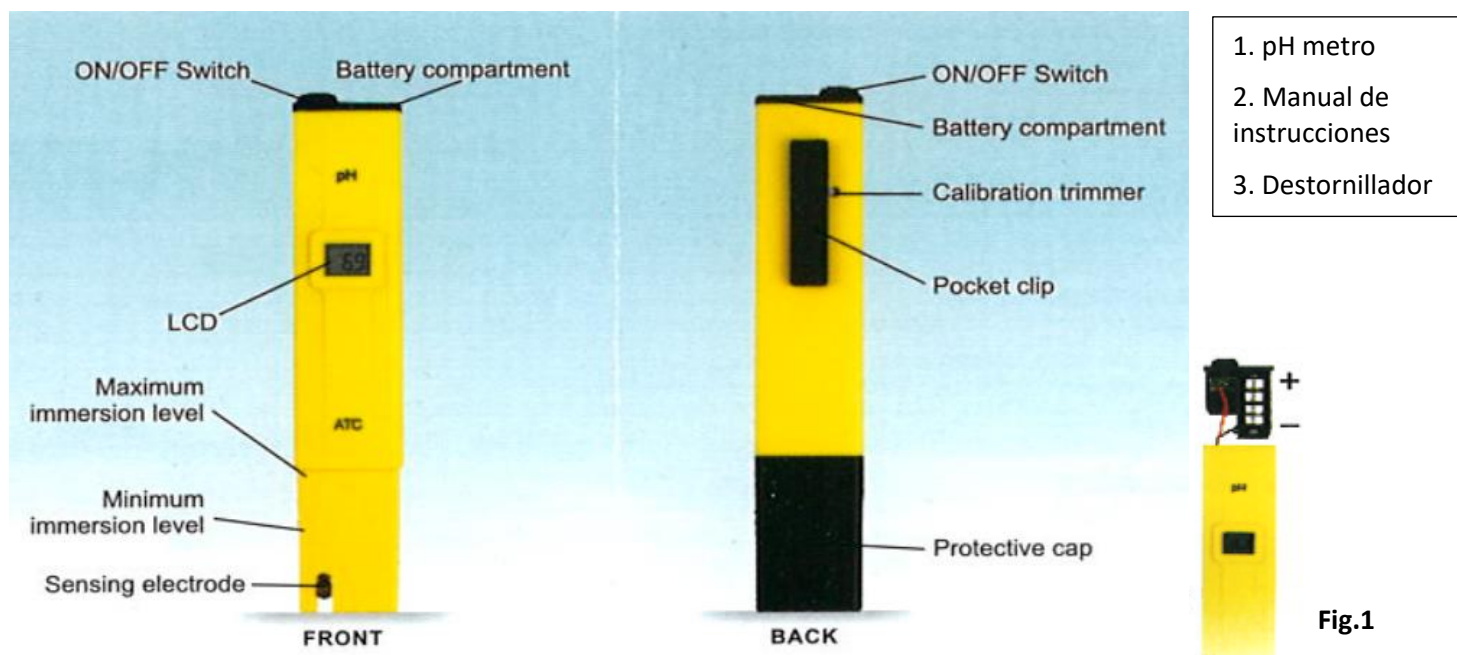
Fig.1

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NOTA: Do not use distilled or deionized water for storage purposes.

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